

1810-B10 and 1810-B20 Bogorov Counting Chamber

Warranty and Parts:

We replace all missing or defective parts free of charge. For additional parts, use part numbers above. We accept MasterCard, Visa, American Express, checks, institutional P.O.'s. All products guaranteed free from defect for 90 days. This guarantee does not include accident, misuse, or normal wear and tear.

Description:

The Wildco[®] Bogorov Chambers are designed to be placed under a binocular dissection microscope in such a way that the groove will be directly under the objective lens. The chamber is constructed of clear acrylic. They are disposable, but you can expect about two to four (2 to 4) weeks of daily use from each one.

Additional Materials

Needed:

- Binocular dissection microscope
- Entomological (curved) forceps (Wildco[®] 7905-T10)
- Entomological probes (Wildco[®] 7905-T25)
- Sample vials (Wildco[®] 7913-B60- 6 oz)
- Disposable pipettes (Wildco[®] 7905-T10)

Warning:

Do not use chemicals such as acetone, ketone and alcohol for cleaning the Bogorov chambers. The use of these chemicals may cause cracking and/or crazing. To remove surface scratches, buff gently with mild abrasive and a soft flannel buffer, then wash as directed.

Rinse preserved samples thoroughly with water before placing them in the Bogorov Chamber. This can be done using a fine sieve and then suspending the samples in water. This will help preserve the chamber and prevent inhalation of the preservative.

How To Use:

1. Place under a low power stereoscope so that the groove is directly under the objective lens. Adjust the stereoscope as necessary to get the best view of the groove.
2. Place a measured sample into the groove using a pipette.
3. Adjust the chamber so the top left-hand end of the groove is in your field of vision.
4. Process your sample by moving the Bogorov Chamber until you reach the other end of the groove.
5. After each use, wash the chamber with mild soap and soft water. Rinse completely with distilled water before drying.

**Order online at
www.wildco.com**

P/N 24-1810

© Wildlife Supply Co. Wildco[®] and Wildlife Supply Co. are registered trademarks of Morris & Lee Inc. All rights reserved.

1810-B10	Bogorov Counting Chamber
1810-B20	Modified Bogorov Counting Chamber
7905-T10	Entomological curved forceps
7905-T25	Entomological probes
7913-B60	Plastic sample vials, wide mouth, 6 oz. capacity
7905-P10	Disposable plastic pipettes